## SEPARATING A MIXTURE BY DISTILLATION

STUDENT BOOK	Chapter 1, page 23	
TOOL BOX	Page 34	
TOOLBOX	raye 34	

## Goal

Separate the constituent substances of a liquid mixture based on their different boiling points.

## **Materials**


- cobalt chloride paper strips
- 50 mL of a liquid mixture
- 125-mL Erlenmeyer flask and two-hole stopper
- 2 or 3 porous stones
- · thermometer
- · hot plate

- condenser tube
- ring stand
- universal clamp
- glass elbow tube and one-hole stopper
- 2 flexible tubings
- 2 100-mL beakers

## **Procedure**



- 1. Dip a cobalt chloride paper strip into the liquid mixture. Record the result.
- 2. Pour the liquid mixture into the flask and add 2 or 3 porous stones.
- 3. Seat the two-hole stopper in the opening of the flask.
- 4. Insert the thermometer into one hole of the stopper so the bulb is submerged completely and not touching the flask.
- 5. Place the flask on the hot plate.
- 6. Secure the condenser tube to the ring stand with the universal clamp.
- 7. Seat the one-hole stopper in the opening at the wide end of the condenser tube and connect the condenser tube to the flask with the glass elbow tube.
- 8. Use flexible tubing to connect the condenser tube water inlet (furthest from liquid mixture) to a cold-water faucet.
- 9. Use flexible tubing to connect the condenser tube water outlet (closest to liquid mixture) to a sink basin.
- 10. Position one beaker under the opening at the narrow end of the condenser tube.
- 111. Open the cold-water faucet. Ensure there is no leak and water flows well into the sink basin.
- 12. Heat the liquid mixture to the boiling point of one constituent substance. Record the boiling point.

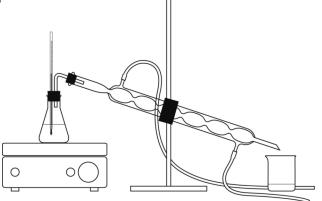

1

Name:	G	oup:	Date:
<ul> <li>13. Continue to heat as long</li> <li>14. Replace the beaker with</li> <li>15. Record the temperature</li> <li>16. Stop heating when evap</li> <li>17. Dip a cobalt chloride page</li> <li>18. Clean up and put away</li> </ul>	the second beaker wat the boiling point. Foration is almost comper strip into each liquidaterials.	when the temperant plete. uid collected. Rec	
<b>Results</b> Record your results in the tab	ble below. Give the ta	ble a title.	
Fitle:			
Characteristic	Liquid mixture	Liquid 1	Liquid 2
<b>1.</b> Were the constituent sub Explain your answer.	-	mixture successfu	ılly separated?
2. According to your results	, what is the nature o	f the liquids colled	cted?
2. According to your results  3. Is it possible to identify the			

5. How could you improve the protocol for this lab?